


MEMORY MONITORING THERMOMETER ALARM INSTRUCTIONS

SPECIFICATIONS

Range:	-58 to 158.0°F / -50 to 70.0 °C
Resolution:	1°
Accuracy:	± 1 °C
Sampling Rate:	Fast 10 seconds or Normal 60 Seconds (user selectable)
Features:	LCD digital thermometer, internal or external display with maximum/minimum memory display and high/low alarm
Battery:	AA Alkaline

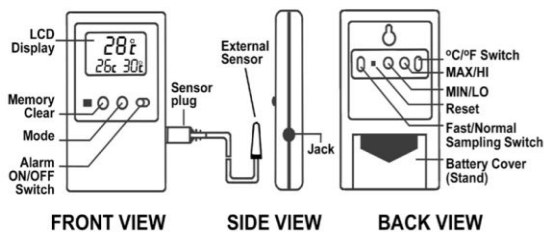

PROBE

The probe is sealed in a miniature bottle (1 x 2-1/2 inches) filled with a patented nontoxic glycol. Solution is GRAS (generally recognised as safe) by the FDA (Food and Drug Administration). It eliminates concerns about incidental contact with food or drinking water. The solution filled bottle simulates the temperature of other stored liquids.

Velcro® and magnetic strip are supplied to mount the bottle to the inside of a refrigerator/freezer and to mount the display unit to the outside. Micro-cable permits refrigerator doors to close on it.

NOTE: Plug probe sensor into the jack on the side of the unit. Press the RESET button once whenever the probe sensor is removed or inserted. Updated ambient or probe temperatures will be displayed. With probe sensor attached the unit simultaneously displays current probe temperature and probe minimum/maximum temperatures.

With the probe sensor removed the unit simultaneously displays current ambient temperature and ambient minimum/maximum temperatures.

RESETTING THE UNIT

Use a pointed object to push the RESET button on the back of the unit. It is necessary to reset the unit whenever a change is made to C/F temperature, FAST/NORMAL sampling rate, or the probe sensor is plugged in or removed.

NOTE: HI/LO temperature alarm settings and MIN/MAX memories will be cleared after resetting the unit.

OPERATION MODES

The unit has two operating modes: NORMAL DISPLAY MODE and ALARM DISPLAY MODE. To change from one mode to another press the MODE button once.

NORMAL DISPLAY MODE – is indicated by the display of current temperature and the MIN (minimum) and MAX (maximum) temperatures. These temperatures relate to external (probe) temperature if the probe is plugged in and to internal (ambient) temperature if the sensor probe is removed.

ALARM DISPLAY MODE – is indicated by the display of current temperature and the HI and LO alarm settings. These temperatures relate to external (probe) temperature if the probe is plugged in and to internal (ambient) temperature if the sensor probe is removed.

RESETTING MIN/MAX MEMORIES

There are two ways to reset the memories:

1. To reset both minimum and maximum memories – press MEMORY CLEAR button.
2. To reset individual minimum and maximum memories – press either the MIN or MAX button (on back of unit) once to reset the respective memory. The other memory will remain unaffected.

NOTE: You must be in the Normal Display Mode to reset the memories.

DISPLAYING °F OR °C

To display the temperature readings in Fahrenheit (°F) or Celsius (°C) slide the switch on the back of the unit to °F or °C.

NOTE: The unit must be reset after changing the temperature unit.

SELECTING THE SAMPLING RATE

To select the normal sampling rate (60 seconds) or fast sampling rate (10 seconds) slide the switch on back of unit to NORMAL or FAST.

NOTE: The unit must be reset after changing the sampling rate. Using the NORMAL sampling rate will extend battery life.

SETTING THE HIGH/LOW ALARM

1. Press the MODE button unit until is in the ALARM DISPLAY MODE indicated by the display of current temperature and HI and LO alarm settings.

2. To set the HIGH temperature alarm limit press the HI button, to advance the display in 1° increments, until the desired high temperature is displayed. Press and hold the HI button to rapidly advance the display in 5° increments.
3. To set LOW temperature alarm limit press the LO button, to advance the display in 1° increments until the desired low temperature is displayed. Press and hold the LO button to rapidly advance the display in 5° increments.

ALARM

The unit will sound the alarm in both NORMAL DISPLAY MODE and ALARM DISPLAY MODE.

Slide the ALARM ON/OFF switch to ON. When the current temperature display rises above the HI set point, or falls below the LO set point, the alarm will sound for 1 minute and issue a 3 second repeater beep every minute thereafter, for up to 12 hours. The alarm will continue to sound even if the current temperature display returns to an “in range” condition.

To temporarily silence the alarm press either the HI or LO button on the back of the unit. The unit is still active and the alarm will sound again if the temperature goes out of range. To permanently silence the alarm, slide the ALARM ON/OFF switch to OFF position.

BENCH STAND

The unit is supplied with a bench stand that is a part of the battery cover. To use the bench stand locate the small rectangular opening at the bottom of the unit. Place your fingernail into the opening and flip the stand up. To close the stand, simply snap it shut.

WALL MOUNTING THE UNIT

Set a screw into the wall at the desired location. Do not set the screw flush to the wall; the head of the screw will need to slip into the receptacle on the back of the unit. Once the screw has been properly set, hang the unit in place by sliding the receptacle on the back of the unit over the head of the screw.

ALL OPERATIONAL DIFFICULTIES

If this thermometer does not function properly for any reason, please replace the battery with a new high quality battery (see “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties.

Replacing the battery with a new fresh battery will solve most difficulties. If the voltage of the battery becomes low °C and °F symbols will flash.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indications the battery must be replaced. Slide the battery cover toward the end of the unit. Remove the exhausted batteries and replace with 2 AA alkaline batteries. Replace the battery cover.

WARRANTY, SERVICE, RECALIBRATION

For warranty, service or recalibration contact:



United Kingdom Office

Woodley Equipment Company Limited
Locomotive House
Locomotive Industrial Estate
Horwich, BL6 5QB
Bolton
United Kingdom
Tel: +44 (0)8456 777001
Fax: +44 (0) 8456 777002



USA Office

Woodley Equipment Company Inc
3718 Northern Boulevard, Suite 512
Long Island City, 11101
New York
USA
Freephone: 1-800-471-9200
Tel: 1-718-606-0516
Fax: 1-718-606-0955

www.woodleyequipment.com

Email: enquiries@woodleyequipment.com

Product Code: WD7TEM001
June 2016