# Vortex Mixer

## SA7 & SA8


Instructions for use

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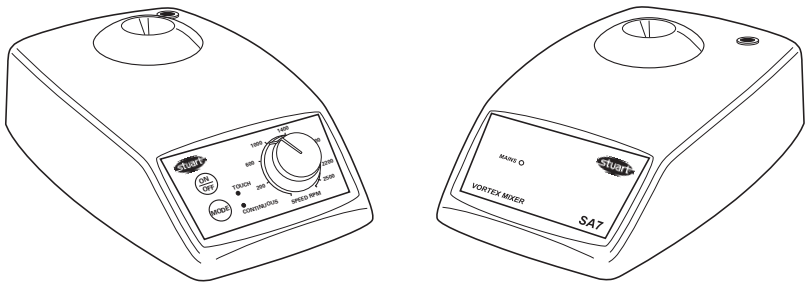

# English

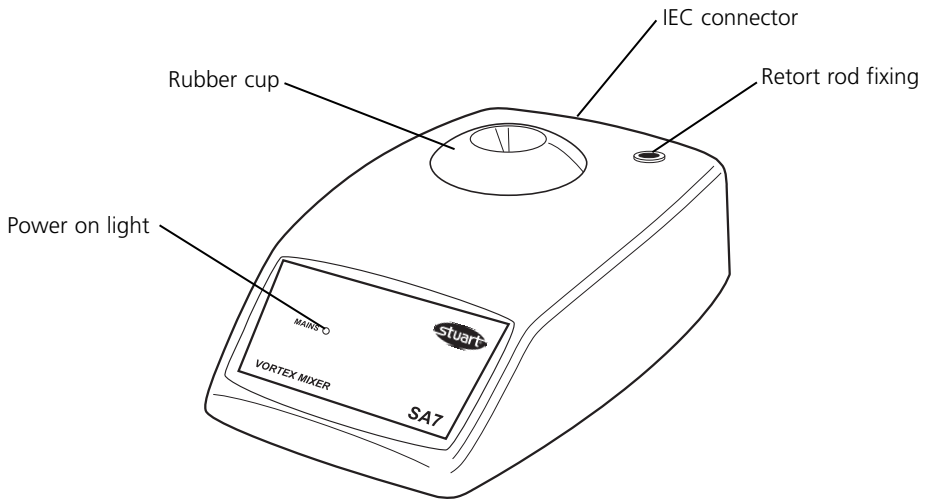**Figure 1: SA7 Fixed Speed model**

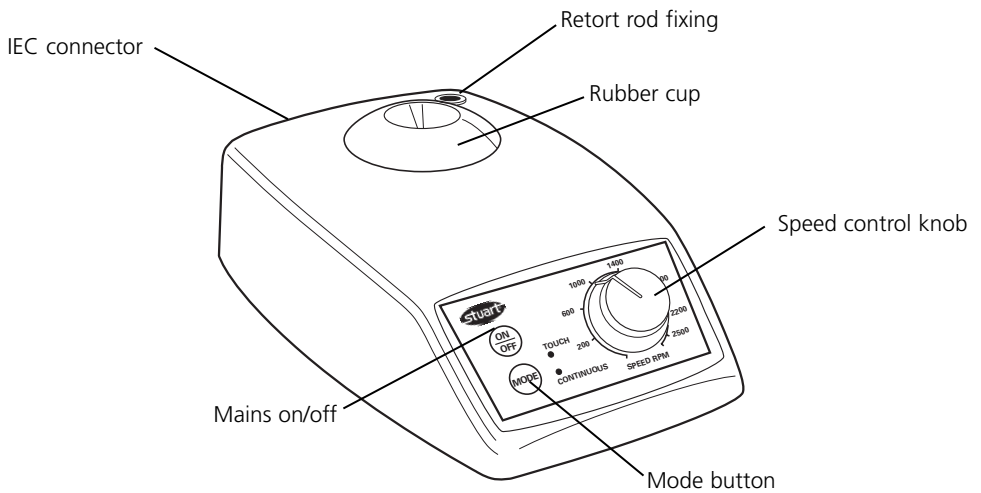**Figure 2: SA8 Variable Speed model**

**Thank you for purchasing this Bibby Scientific product. To get the best performance from the equipment, and for your personal safety, please read these instructions carefully before use.**

This equipment is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuations not exceeding 10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all round of 200mm from walls or other items

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

## Electrical Installation

**THIS EQUIPMENT MUST BE EARTHED**

**Before connection please ensure that the line supply corresponds to that stated on the rating label.**

### Power requirements:

Model	Wattage
SA7	20W
SA8	20W

There is an IEC socket at the rear of the instrument for connection to the mains supply, (see figure 2). The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a U.K. 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable, take the lead with the U.K. plug and replace the plug with a

suitable alternative. See the enclosed instruction sheet for advice on how to carry out this procedure.

Should the mains lead require replacement a cable of 1mm<sup>2</sup> of harmonised code H05W-F connected to an IEC 320 plug should be used.

### **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN**

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

### **Safety Advice Before Use**

- ❖ Never move or carry the unit when in use or connected to the mains electricity supply.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In the case of mains interrupting, a fault or mechanical failure, the unit will NOT continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

## Operation

Position the apparatus on a firm level surface, ensuring that the IEC connector is in easy reach.

### **SA7 Fixed speed version (see Fig 1)**

Connect the unit to the mains. Green neon lights. To mix tube, press base firmly down onto the rubber cap, the unit will start. Once the contents of tube are mixed, lift clear, the unit will stop until another tube is pressed onto the cup.

### **SA8 Variable speed version (see fig 2)**

Ensure that the speed control knob is fully to the left. Connect the unit to the mains.

Turn ON by pressing the <On / Off> button once (a green neon lights to show the unit is in the touch mode).

To mix tube, press base firmly down onto the rubber cap. Once the contents of tube are mixed, lift clear. Select the optimum mixing speed by turning the control knob to the right.

To select continuous mode, press the <mode> button (the light will change to show continuous mode). The unit will now remain on until either the <mode> button is pressed to select touch mode or the unit is switched OFF.

### Accessory heads

**N.B. Accessory heads are for use with model SA8 only. Accessory heads should always be used in the continuous mode.**

To change the head, remove the rubber cup by pulling upwards and store in a safe place. Insert the accessory head ensuring that the two slots in the head are in line with the lugs in the vortex mixer top.

The microtitre plate holder (SA8/1) should not be used at speeds above 800rpm.

## Maintenance & Servicing

 **WARNING:** Ensure the unit is disconnected from the mains electricity supply before attempting maintenance or servicing.

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

**Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.**

The following spares and accessories are available from your laboratory dealer.

Description	Catalogue Number
Microtitre plate holder	SA8/1
Spare rubber cup	SA7013
Retort rod, 300mm	SR2

For a comprehensive list of parts required by service engineers conducting internal repairs and a service manual, please contact the Technical Service Department of Bibby Scientific Ltd. quoting both the model and serial number.

Only spare parts supplied or specified by Bibby Scientific Ltd. or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

If in any doubt, please contact the Technical Service Department of Bibby Scientific Ltd. or the point of sale.

Bibby Scientific Ltd.  
Stone, Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)844 936 0234  
Fax: +44 (0)844 936 0235  
e-mail [info@bibby-scientific.com](mailto:info@bibby-scientific.com)  
[www.bibby-scientific.com](http://www.bibby-scientific.com)


## Warranty

Bibby Scientific Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of three (3) years. In the event of a justified claim Bibby Scientific will replace any defective component or replace the unit free of charge.


This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non approved parts or repair by unauthorised personnel.

# Technical Specification

	<b>SA7</b>	<b>SA8</b>
Speed, rpm	2,500	250-2,500
Orbit, mm	4.2	4.2
Continuous mode	No	Yes
Accessory heads	No	Yes
Maximum load	1 Kg	
Overall dimension	78 x 135 x 215mm	
Net weight	3.2 kg	
Electrical supply	90 - 240V, 20W	

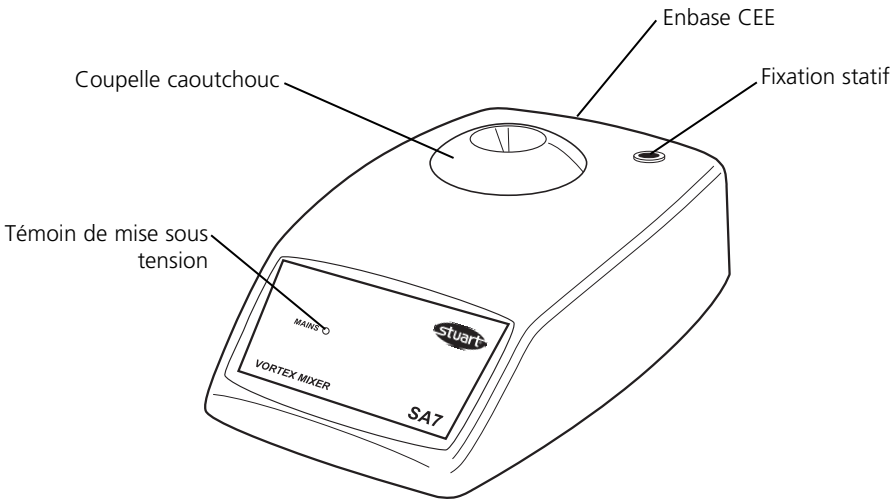**Fig. 1: Version vitesse constante**

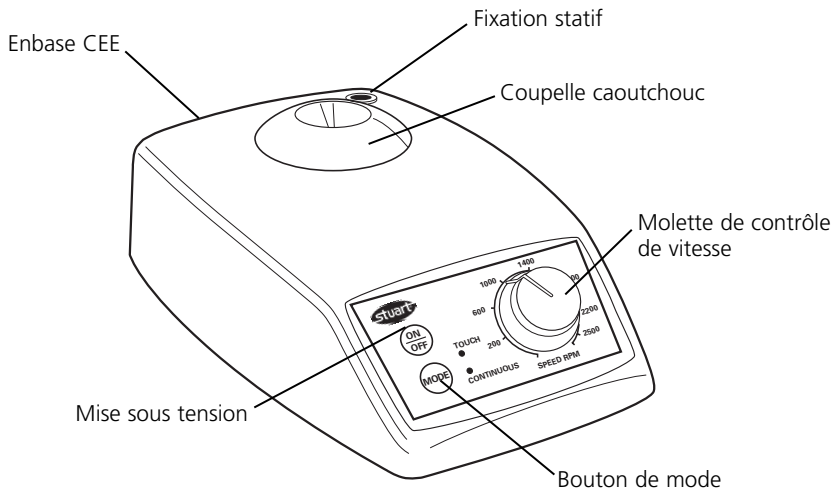**Fig. 2: Version vitesse variable**

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# Notes

These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Scientific Ltd.

## *Declaration of Conformity*

---

### **Vortex Mixer**

#### **Model SA7 & SA8**

These products comply with the requirements of the EU Directives listed below:

**89/336/EEC** Electromagnetic Compatibility Directive, amended by 92/31/EEC and 93/68/EEC.

**73/23/EEC** Low Voltage Directive amended by 93/68/EEC

Compliance with the requirements of these Directives is claimed by meeting the following standards:


EN 61326: 1997 + Amendments A1: 1998, EN 61000-4-2: 1995, EN 61000-4-4: 1995, EN 61000-4-5: 1995, EN 61000-4-6: 1996, EN 61000-4-9: 1993, EN 61000-4-11: 1994, EN 61000-3-2: 1995, EN 61000-3-3: 1995, EN 61010-1: 1993 + Amendments A1: 1992 & A2: 1995, EN 61010-2-51: 1995.

#### Compliance Certificates and Full Reports.

Ref: RETS0287/31 and RETS0287/32

From "Epsilon Technical Services" an Independent Accredited Test House Showing Compliance to the above Standards are available on request.

**CE mark affixed '01**

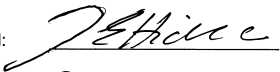Signed:  (Mr D. E. Hicks)

Date: 2nd December 2005

Authority: Technical/Development Manager

for

Barloworld Scientific Ltd - Stone - Staffs - ST15 0SA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

# INSPECTION REPORT

**MODEL** \_\_\_\_\_


## ELECTRICAL SAFETY

- 1. Earth continuity
- 2. Insulation
- 3. Flash test

## FUNCTIONAL

- 1. Indicators
- 2. Vortex action
- 3. Visual acceptance

**QUALITY CONTROL INSPECTOR**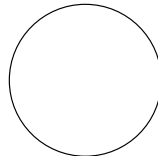

### **Bibby Scientific France SAS**

ZI du Rocher Vert - BP 79  
77793 Nemours Cedex  
France  
Tel: +33 1 64 45 13 13  
Fax: +33 1 64 45 13 00  
e-mail: bsf@bibby-scientific.fr  
www.bibby-scientific.com

### **Bibby Scientific Italia Srl**

Via Alcide de Gasperi 56  
20077 Riozzo di Cerro al Lambro  
Milano Italia  
Tel: +39 (0)2 98230679  
Fax: +39 (0)2 98230211  
e-mail: marketing@bibby-scientific.it  
www.bibby-scientific.it

### **Bibby Scientific Ltd**

Beacon Road Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)844 936 0234  
Fax: +44 (0)844 936 0235  
email: info@bibby-scientific.com  
www.bibby-scientific.com

### **Bibby Scientific US Ltd**

3 Terri Lane Suite 10  
Burlington NJ 08016  
USA  
Tel: 800-225-9243  
Fax: 609-589-2571  
www.bibby-scientific.com

### **Afora S.A.**

Calle Aribau 240  
08006 Barcelona Spain  
Tel: +343 93-306 98 00  
Fax: +343 93-306 98 23  
e-mail: marketing@afora.com  
www.afora.com